


## Prepare your slide

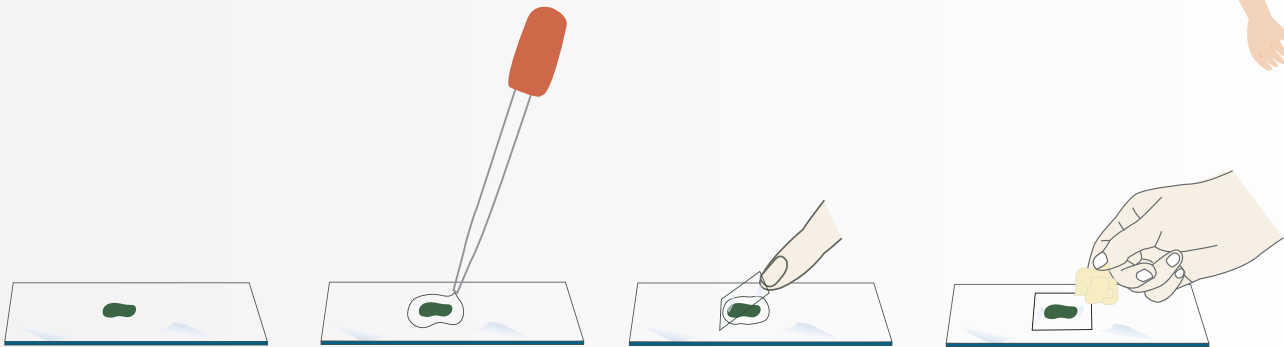

### Examine your water sample

You can't wait to examine the water from your garden or a park? Then take a water sample and let us prepare for your microscope. With it, you can take a close look at what's swimming around.

To do this, take a microscope slide and put one or two drops of your sample on it with a pipette (eye dropper). Then put the thin cover glass over it. If necessary, press down the cover glass lightly. You can gently soak up the excess water with paper towel on one side.

Do you want to colour your finished sample? Use a drop from a normal ink cartridge, e.g. from your fountain pen. Put the ink on the left or right side edge of the cover glass. Hold a blotting paper or filter paper on the opposite edge. You will see: The ink will be drawn through your entire sample and stain it. Now you are ready to examine your first specimen - let's go!



Find us on:



[www.bresser-junior.com](http://www.bresser-junior.com)